• Gram positive spherical or ovoid cells, 0.7-0.9 µm in diameter in pairs to long chains. Final pH range in glucose broth is 4.0-4.5. The preferential fermentation of the disaccharides, sucrose and lactose, may result in a lower pH value as compared to glucose fermentation. Acid is produced from glucose, fructose, lactose and sucrose; no acid from trehalose, maltose, inulin, glycerol, mannitol, sorbitol or salicin and rarely from raffinose, xylose or arabinose. Optimum temperature is between 40°C and 45°C. Growth occurs at 50°C but not at 53°C. No growth at temperatures below 20°C. Heat tolerance: survives 65°C for 30 min. Source: milk and milk products such as cheese and yoghurt. Often used as a starter culture for these products. This species is easily recognized by its thermal tolerance; unable to ferment maltose and unable to grow in media containing ≥2.0% sodium chloride.

Homemade Yoghurt (a recipe for your own houshold as well)	1. Inoculum is a tenth of the volume (with one 2 dl cup of white yoghurt you can inoculate 2 l of milk).
	2. Either boil the milk shortly or better keep it at 85°C for 30 min. Then let it cool to 45°C.
	3. Add the inoculum while stirring the milk.
	4. Fill milk-yoghurt mixture into clean glasses, cover with aluminium foil.
	5. Heat water to 45°C in a big pan, take pan from the hot plate, put glasses inside and cover the pan with a heat insulation (a woolen blanket for instance), or incubate the glasses in a 45°C preheated oven.
	6. Leave at 45°C over night, then put into the refrigerator.
	Remark: The initial heating to 85°C during 30 minutes denatures some of the milk proteins which afterwards leads to a yoghurt of a more stable and finer consistency. At 45°C, the yoghurt formation is finished after about 4 hours, longer incubation times will lead to a more solid and more acidic tasting yoghurt.
Material	Each group gets: - 6 clean, sterile (autoclaved) glass beakers, 100 ml, covered with aluminium foil
	- 120 ml of milk - 1 teaspoon
	Available:
	- plain yoghurt bought in a food store
	- tripod, wire mesh, Bunsen burner
	 paper towels 4°C, 37°C and 45°C incubators
	- pH paper
Laboratory Rules &	No microbial risks involved.
Precautions	Keep the inoculum uncontaminated. Watch milk carefully while heating it.
Experiences gained	Taste a little bit of microbiology!
Timing	30 minutes
Reporting	Explain the purpose of the 8 steps outlined in the Experimental Protocol. Take notes on the experiment in your laboratory journal and report the results in class.